

Pro-GA <Cell-permeable GGCT inhibitor>

Catalog NO. FDV-0019

Research use only, not for human or animal therapeutic or diagnostic use.

Product Background

γ -Glutamylcyclotransferase (GGCT) catalyzes the formation of 5-oxoproline and free amino acids from γ -glutamyl dipeptides for maintaining glutathione homeostasis. Although the enzymatic activity of GGCT was firstly reported in 1956, the gene of GGCT had not been identified for many years. In 2008, a biochemical study revealed C7orf24, a functionally unknown and highly expressed protein in bladder cancer cells, shows GGCT activity (ref.1, 2). Many studies show GGCT is highly expressed in various cancer cells as compared with normal tissues. Overexpression of GGCT promotes cell proliferation in NIH-3T3 with low-level GGCT expression and depletion of GGCT mRNA by RNA interference inhibits cell-growth in several types of cancer cell lines which show high expression of GGCT. A potent inhibitor of GGCT like RNAi methods will be powerful tool to investigate the physiological function of GGCT. *N*-Glutaryl-*L*-alanine (GA) is one of the potent inhibitor for GGCT activity, but GA has little membrane-permeability. Pro-GA is a novel diester-type cell-permeable GGCT inhibitor pro-drug and shows the inhibitory effects for proliferation of multiple malignant cancer cells *in vitro* and tumor growth *in vivo* (Fig.2, ref.3).

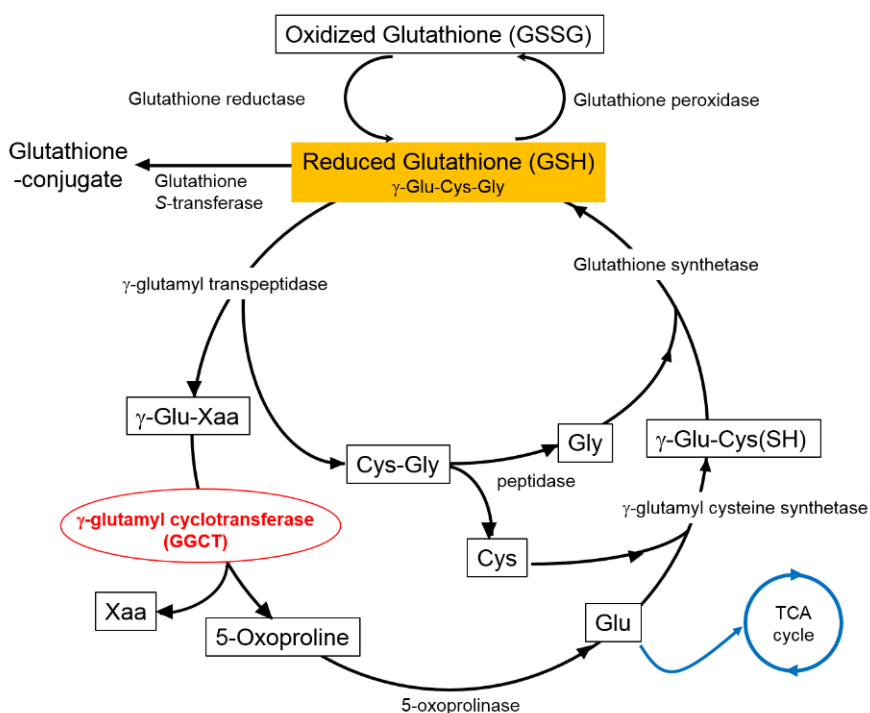


Figure 1. Overview of glutathione metabolic cycle

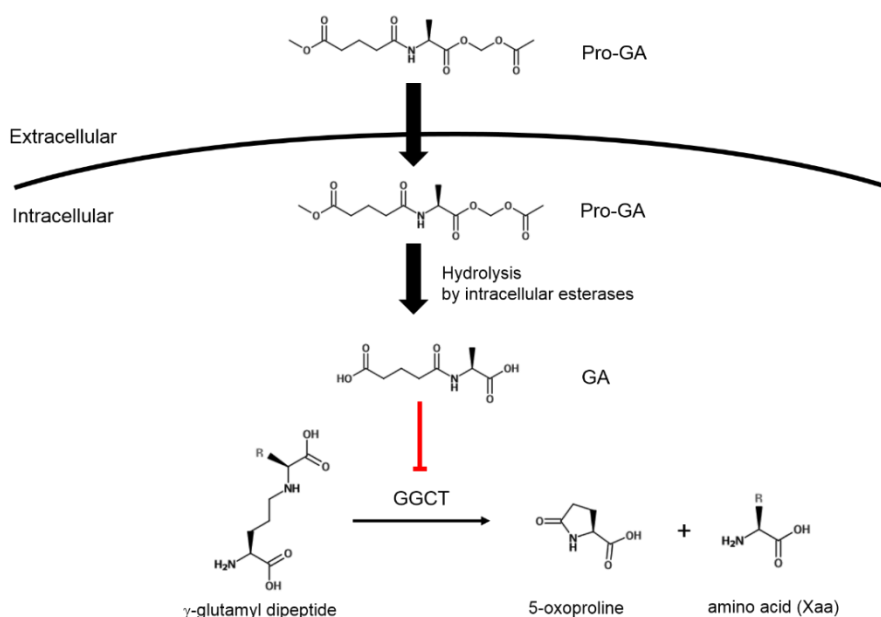


Figure 2. Principle of pro-drug type inhibitor Pro-GA for GGCT

Description

Catalog Number: FDV-0019
 Size : 2 mg
 Formulation : C₁₂H₁₉O₇N
 Molecular weight : 289.28 g/mol
 Solubility : Soluble in DMSO

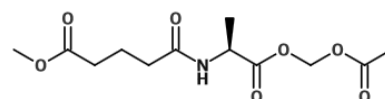


Figure 3. Chemical structure of Pro-GA

Reconstitution and Storage

Reconstitution : Stock solution recommended concentration 0.2-1 M in 100% DMSO.
 Storage (solution) :
 Store powder at -20°C.
 After reconstitution in DMSO, aliquot and store at -20 °C. Avoid repeated freeze-thaw cycles.

How to use

General procedure of cell-based inhibition of GGCT by Pro-GA

<Before Use>

pH of assay medium is very important to avoid non-enzymatic hydrolysis of Pro-GA. Because both acidic and basic conditions can induce hydrolysis of Pro-GA, pH should be neutral. As the assay medium, DMEM containing 25 mM HEPES (DMEM/HEPES) is strongly recommended.

1. Prepare 100 μM of Pro-GA in DMEM/HEPES just before use.

Note: Pro-GA is not stable in DMEM/HEPES for long time.

Please use Pro-GA in DMEM/HEPES within 5 min.


2. Exchange culture medium to the Pro-GA in DMEM/HEPES and culture cells for 1-120 hours.

Note: Pro-GA is immediately penetrated into cells and converted to GA inside the cells by esterases. Regarding detail information, please find “GA-release efficiency of Pro-GA in cells” below.


Application data

GGCT inhibition activity of parent compound GA *in vitro*

GGCT activities in cell lysates of NIH3T3 or GGCT-overexpressed NIH3T3 were measured by fluorescent substrate LISA-101. 100 μ M GA strongly inhibited endogenously expressed GGCT activity.

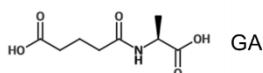


Note: This data is only for reference of an inhibition activity of GA *in vitro*. Pro-GA could not be used *in vitro* assay such as enzymatic activity in the cell lysate.

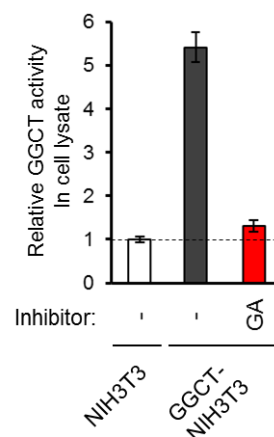


Pro-GA inhibits the proliferation of cancer cells

Indicated normal cells and cancer cells were cultured in serum starve condition with DMSO (as control), GA and Pro-GA for the indicated number of days. Cell proliferation was calculated by the WST cell viability assay. Pro-GA specifically inhibited proliferation of cancer cells.

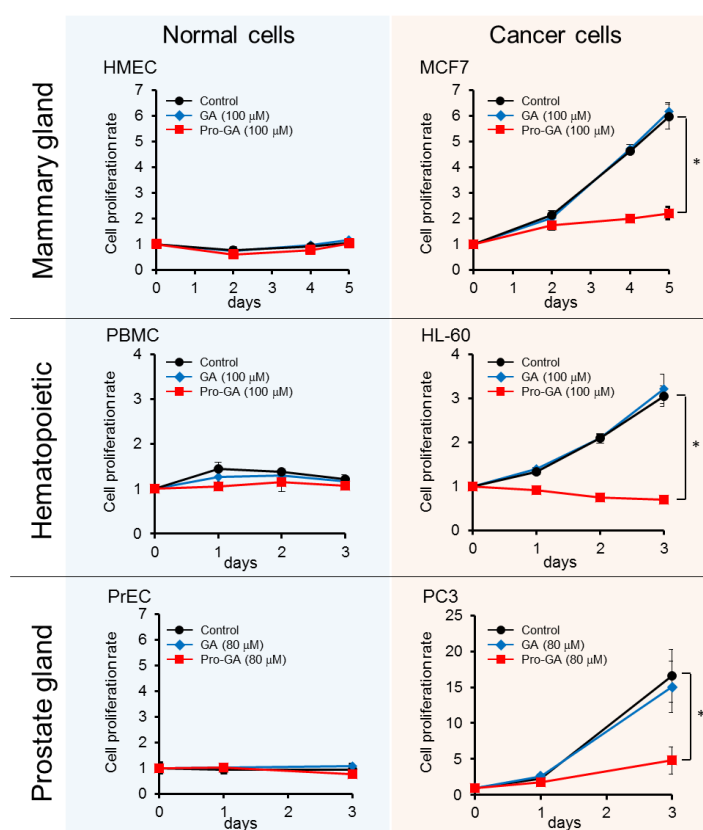


Pro-GA inhibits tumor growth in xenograft mouse model

PC3 tumor cells were implanted into CB17 SCID mice and DMSO (as control) or Pro-GA was intraperitoneally administrated twice a week (n=5 per group). Administration of Pro-GA significantly inhibited tumor growth *in vivo*.

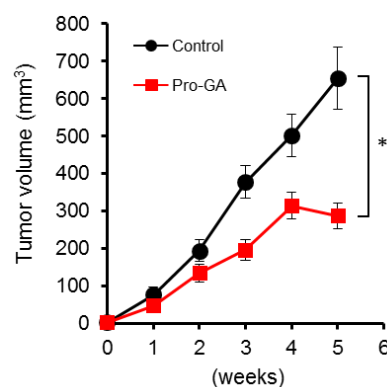


Stability of Pro-GA in culture medium at room temperature

Stabilities of fluorescent NBD-conjugated Pro-GA in PBS or 10% FBS/DMEM were monitored by HPLC analysis. While conversion of Pro-GA to GA were slowly in PBS ($t_{1/2}$ >5 hours), Pro-GA was quickly converted into GA in 10% FBS/DMEM by esterase activity in FBS ($t_{1/2}$ ~50 min). Please prepare Pro-GA solution in assay medium just before use.

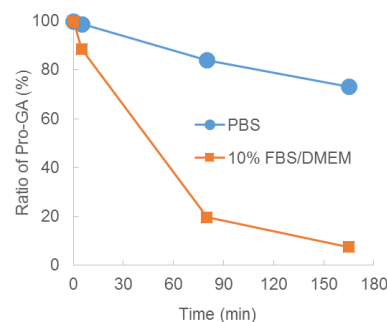

GA-release efficiency of Pro-GA in cells

MCF7 cells were treated with 100 μ M of fluorescent NBD-conjugated Pro-GA for 0-60 min. After wash cells, cell lysates were prepared and analyzed by HPLC to evaluate GA amount released from Pro-GA. NBD-conjugated Pro-GA was immediately incorporated into cells within a few minutes and converted to GA in the cells.

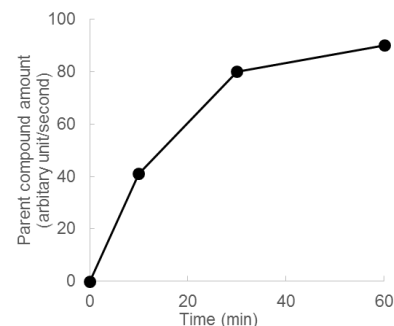

Reference

1. Oakley *et al.*, *J. Biol. Chem.* **283**, 22031-22042 (2008) The identification and structural characterization of C7orf24 as γ -glutamyl cyclotransferase an essential enzyme in the γ -glutamyl cycle
2. Kageyama *et al.*, *Proteomics Clin. Appl.*, **1**, 192-199 (2007) A novel tumor - related protein, C7orf24, identified by proteome differential display of bladder urothelial carcinoma
3. Ii *et al.*, *ChemMedChem.*, **13**, 155-163 (2018) A Novel Prodrug of a γ -Glutamylcyclotransferase Inhibitor Suppresses Cancer Cell Proliferation in vitro and Inhibits Tumor Growth in a Xenograft Mouse Model of Prostate Cancer

Disclaimer/免責事項

This product has been commercialized by Funakoshi Co., Ltd. based on the results of academic research, and the advertisement text, figures and manuals (hereinafter "Product information") have been prepared based on published research reports on March, 2018. The academic interpretation at the time of creation of the Product Information may change in accordance with future developments in the relevant research field and expansion of various scientific findings, and the latest version and certainty of the Product Information are not guaranteed. The specifications of this product and the Product Information are subject to change without notice. Please contact us for the latest information.

本商品は学術研究成果を基にフナコシ株式会社が商品化したもので、2018年3月時点における公開研究報告を基に広告文章およびマニュアル(以下、商品資料)を作成しています。今後の当該研究分野の発展および各種学術知見の拡大にともない、商品資料作成時の学術的解釈が変更になる可能性があり、最新性・確実性を保証するものではありません。また、本商品の仕様および商品資料を予告なく変更する場合がございます。最新の情報に関しましては、弊社までご確認いただきますようお願い申し上げます。

E-mail Newsletter Sign Up

Japanese English

 **funakoshi**
FRONTIERS IN LIFE SCIENCE
URL: <http://funakoshi.co.jp>
9-7 Hongo 2-Chome, Bunkyo-ku, Tokyo 113-0033