

LentiBlast Premium

INSTRUCTION MANUAL

Lentivirus Transduction Enhancer



OZ BIOSCIENCES
The art of delivery systems



Instruction Manual

Description

LentiBlast Premium The Lentiviral Transduction Enhancer

LentiBlast Premium is a chemical lentiviral transduction enhancer, very effective for enhancing viral driven genetic modification. LentiBlast Premium is non-cytotoxic and improves transduction in a wide range of cell types from CD34+ hematopoietic Stem cells to primary cells or cell lines such as HT1080¹ or SH-SY5². Moreover, its formulation ensures total compatibility and performance for enhancing infection and transduction in CD4+ and CD8+ T lymphocytes^{3,4} and is perfect for CAR-T therapy⁵. The efficiency of LentiBlast Premium is such that the viral MOI (number of viral particles per cell) can be reduced and therefore fewer viruses can be used.

Catalog Number	Description	Volume (µL)	Number of transductions in a 24-well plate (average)
LBPX500	LentiBlast Premium 100 transductions	500 µL	100
LBPX1500	LentiBlast Premium 300 transductions	1500 µL	300

For any technical question, contact us at tech@ozbiosciences.com

1. Technology

1.1. Description

LentiBlast Premium is the ideal reagent to enhance lentiviral infection and transduction in any type of cells: adherent or in suspension, primary or cell lines. Its patented chemical composition allows to simultaneously neutralize electrostatic repulsions between membrane and viral particles and to enhance viral fusion with cell membrane. Due to a favorable “membrane permeable effect” limiting the transmembrane potential changes, LentiBlast Premium is non-toxic.

NOTE: LentiBlast Premium can also be used with other enhancers and for sequential transductions.

1.2. Storage and shipping condition

Storage: Upon reception and for long-term use, store the LentiBlast Premium transfection reagent at -20°C

Stability: 1 year

Shipping condition: The reagent is shipped at RT

¹ Arnoult N., Nature. 2017 (549), 548–552.doi:10.1038/nature24023

² Araújo-Vilar D., Eur J Hum Genet. 2018 Jan 24. doi: 10.1038/s41431-017-0052-8

³ Epstein AL., Patent, 2016

⁴ Benati D., J Clin Invest. 2016 Jun 1;126(6):2093-108.

⁵ Zheng L., Int J Mol Sci. 2017 Dec 20;18(12). pii: E2773

2. Applications and Protocols

2.1. General Considerations

- Allow reagent to reach room temperature before starting
- Do not change the transduction conditions already settled: simply add the LentiBlast Premium to your already established infection/transduction protocol
- If transduction conditions are unknown, we recommend beginning with a MOI of 2
- We suggest to first test 3 dilutions of LentiBlast Premium (1:1000, 1:100 and 1:50)

2.2. Cells preparation

Cell culture prior to transduction: the day before transduction prepare the cells according to the table below.

Effects of lentiviral transduction are generally observed after 48 to 96 h incubation. Cells should be 20-50 % confluent at the time of transduction (see the suggested cell number in the Table 1).

Tissue culture dish format	Surface area per well ¹	Cell number	Recommended volume for culture	Recommended volumes for viral preparation
96-well plate	0.3 cm ²	3 - 8 x 1.10 ³	100 µL	50 µL
24-well plate	2 cm ²	2 – 4 x 1.10 ⁴	500 µL	50 µL
12-well plate	4 cm ²	4 – 8 x 1.10 ⁴	1 mL	100 µL
6-well plate	9 cm ²	1 – 2 x 1.10 ⁵	2 mL	200 µL


¹ Surfaces area may vary depending on the manufacturer.

Table 1: Recommended cell number and volume for viral/LentiBlast Premium complexes preparation

2.3. Standard Protocol

Use the following protocol to find the ideal conditions for LentiBlast Premium in 24-well plate. Adjust volumes to different well formats according to Table 1. If the lentiviral transduction/infection conditions are unknown, we recommend starting with a MOI of 2 using a lentiviral vector encoding for a fluorescent protein.


1. VIRUS PREPARATION

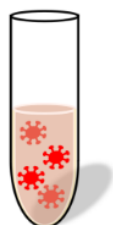

Dilute virus into culture medium sufficient for 4 samples (50 μ L each).

MOI 2 is recommended in case of unknown lentiviral transduction conditions.

2. DISPATCH EQUAL VOLUME OF VIRAL SUSPENSION INTO 4 TUBES

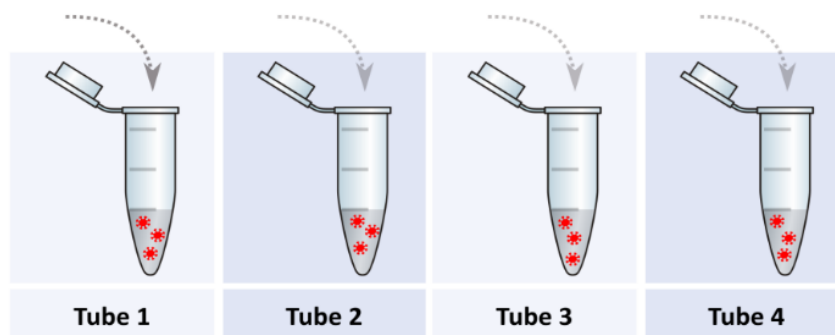

3. ADD LENTIBLAST TO EACH TUBE

	Tube 1	Tube 2	Tube 3	Tube 4
LentiBlast	-	0.5 μ L	5 μ L	10 μ L

4. MIX VIALS BY INVERTING

Do not vortex or centrifuge

5. ADD VIRUS +/- LENTIBLAST


Incubate the cells 24 h under standard culture conditions

6. Optional: 24 h MEDIUM EXCHANGE

Remove medium from the cells
add pre-warmed culture medium

7. INCUBATE CELLS 24 TO 96 h.

Incubate the cells under standard culture conditions
We recommend performing assay from 24 to 96 h.

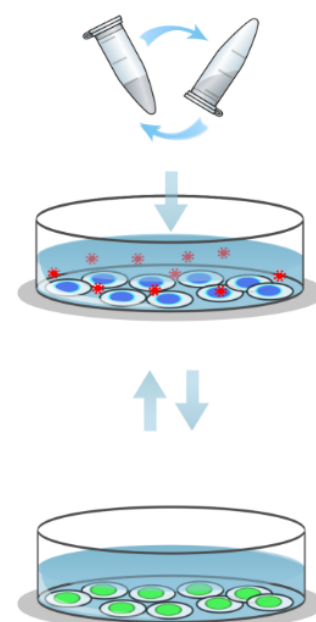

2.4. Optimization protocol

Depending on cell type (primary/cell lines, adherent/suspension, ...), virus serotype, gene of interest, culture conditions, readout..., transduction or infection efficiencies may vary. It may be necessary to optimize the conditions by (1) varying LentiBlast Premium amounts, (2) finely tuning transduction parameters or (3) centrifuging the cells.

2.4.1. Optimizing LentiBlast Premium volume

Using a fixed MOI, vary LentiBlast Premium volumes according to the Table 2 below:

Tissue culture dish format	1:1000	1:500	1:250	1:100	1:50	1:25
96-well plate	0.15 µL	0.3 µL	0.6 µL	1.5 µL	3 µL	6 µL
24-well plate	0.5 µL	1 µL	2 µL	5 µL	10 µL	20 µL
12-well plate	1 µL	2 µL	4 µL	10 µL	20 µL	40 µL
6-well plate	2 µL	4 µL	8 µL	20 µL	40 µL	80 µL

Table 2: Recommended dilutions and volumes of LentiBlast Premium depending on tissue culture dish format

2.4.2. Finely tuning transduction parameters

We have observed that replacing culture medium at the time of transduction with complete medium containing viral particles/LentiBlast Premium could improve the overall efficiency.

- 1) Seed the cells as previously described
- 2) Prepare a viral suspension in complete culture medium; volume should correspond to cell culture volume (refer to table 1 for recommended volumes)
- 3) Add LentiBlast Premium to the viral suspension
- 4) Mix vial by inverting
- 5) Replace cell culture medium with viral suspension/LentiBlast Premium solution
- 6) Optionally perform a medium change after 24H, and incubate 24H to 96H

2.4.3. Raising efficiency using Centrifugation step

For hard-to-transduced cells, a centrifugation step can be added to the standard protocol. Cells are prepared the day of transduction, counted, pelleted and suspended in Lentivirus/LentiBlast Premium mix before centrifugation.

NOTE: Centrifugation may negatively influence cell viability.


- 1) Detach cells and seed them into 4 wells. Refer to Table 1 for suggested cell density
- 2) Follow steps 1 to 4 of the standard protocol and add lentivirus/LentiBlast Premium mixes to cells
- 3) Centrifuge the plate 900 rpm for 90 min
- 4) Incubate cells overnight and proceed to steps 6 and 7 of the standard protocol

3. Troubleshooting

Problems	Comments and Suggestions
Low transduction efficiency	<p>1- MOI used is low. Raise the lentiviral particles amount: up to MOI 100 can be used.</p> <p>2- Infectious viral titer is low. Check viral titer via transduction of HEK-293T cells with serial dilutions of lentivirus.</p> <p>3- Cells are difficult to transduce. Use the centrifugation protocol (refer to paragraph 2.4.3).</p> <p>4- No effect of LentiBlast Premium is observed. Concentrations of LentiBlast Premium is too low or not optimized (refer to paragraph 2.4.1).</p>
Low viability	<p>1- MOI is too high. Lower the lentiviral dose</p> <p>2- LentiBlast Premium concentration is too high. Decrease concentration of LentiBlast Premium, refer to Table 2.</p> <p>3- Cells are sensitive to lentiviral treatment. Perform a medium change after 4 to 6 h or directly after the centrifugation process.</p>

4. Related products

APPLICATIONS

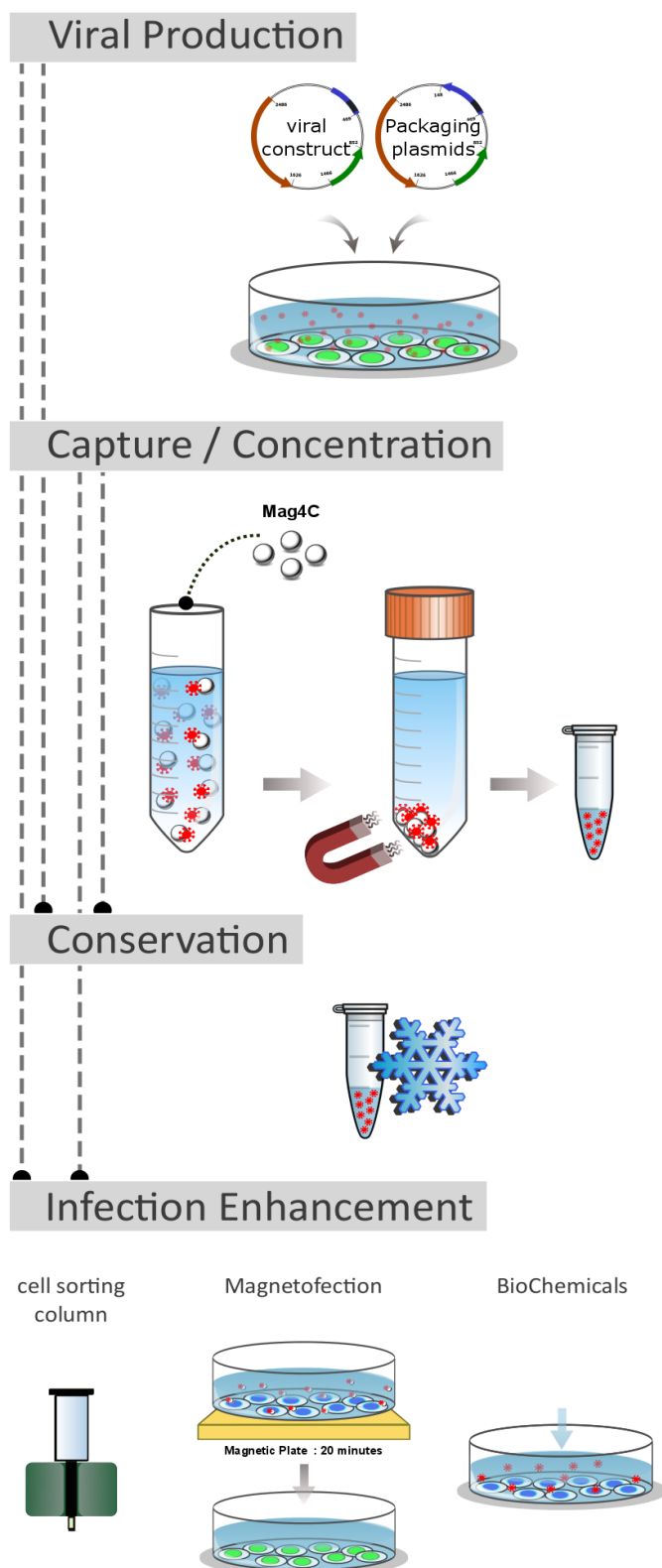

RECOMMENDED PRODUCTS

Calcium Phosphate Transfection Kit

**DreamFect Gold
Helix-IN Transfection Reagent**

High transfection efficiency
High compaction level
Biodegradable

**Mag4C Kits
Mag4C-Lv
Mag4C-Ad**

Concentrate virus
Raise viral dose
Isolate virus from inhibitor
Conserve the viral particles

**Mag4C
conservation buffer**

Adenoviral particles
Retro/lentiviral particles

Magnetic Particles

**Viro-MICST
ViroMag
AdenoMag**

Raise transduction efficiency
Synchronize infection

BioChemicals

**LentiBlast
AdenoBlast**

Enhance transduction
Non toxic
Low MOI

Our technical support will be pleased to answer any of your request and to assist you in your experiments. Do not hesitate to contact us for all complementary information and remember to visit our website in order to stay inform on our latest breakthrough technologies and update:

<http://www.ozbiosciences.com>

Purchaser Notification

Limited License

The purchase price paid for LentiBlast Premium by end users grants them a non-transferable, non-exclusive license to use the kit and/or its separate and included components (as listed in the section 1, Kit Contents). These reagents are intended **for internal research only** by the buyer. Such use is limited to the use in the product manual. Furthermore, research only use means that this kit and all of its contents are excluded, without limitation, from resale, repackaging, or use for the making or selling of any commercial product or service without the written approval of OZ Biosciences.

Purchasers may terminate this License at any time by returning all LentiBlast Premium material and documentation to OZ Biosciences, or by destroying all LentiBlast Premium components. Purchasers are advised to contact OZ Biosciences with the notification that a LentiBlast Premium is being returned in order to obtain a refund and/or to expressly terminate a research only license granted through the purchase of the kit(s).

This document covers in full the terms of the LentiBlast Premium research only license, and does not grant any other express or implied license. The laws of the French Government shall govern the interpretation and enforcement of the terms of this License.

Product Use Limitations

The LentiBlast Premium is developed, designed, intended, and sold for research use only. It is not to be used for human diagnostic or included/used in any drug intended for human use. All care and attention should be exercised in the handling of the kit components by following appropriate research lab practices.

For more information, or for any comments on the terms and conditions of this License, please contact:

Director of Business Development OZ Biosciences SAS

Parc Scientifique et Technologique de Luminy
Bâtiment Grand Luminy technopole
Case 922 zone entreprises
13288 Marseille Cedex 9, France
Ph: +33 (0)4.86.94.85.16
Fax: +33 (0)4.86.94.85.15
E-mail: contact@ozbiosciences.com

EUROPE & ASIA OZ Biosciences SAS

163 avenue de Luminy
Case 922, zone entreprise
13288 Marseille cedex 09
France
Ph: +33 (0) 486 948 516
Fax: +33 (0) 486 948 515
contact@ozbiosciences.com
order@ozbiosciences.com
tech@ozbiosciences.com



USA & CANADA OZ Biosciences INC

4901 Morena Blvd
Suite 901
San Diego CA 92117
USA
Ph: + 1-858-246-7840
Fax: + 1-855-631-0626
contactUSA@ozbiosciences.com
orderUSA@ozbiosciences.com
techUSA@ozbiosciences.com