

# WST-8 Cell Proliferation Kit

## INSTRUCTION MANUAL

### Assays kits



### *WST-8 Cell Proliferation Kit*

The WST-8 Cell Proliferation Kit is a ready-to-use assay kit for determination of viable cell number and for studying induction or inhibition of cell proliferation *in vitro*.


	Content	Catalog Number	Number of assays (96-well plate)
<b>WS1000</b>	10 mL of WST-8 Solution	WS1000	1000

For any technical questions, contact us at [tech@ozbiosciences.com](mailto:tech@ozbiosciences.com)

# 1. Technology

## 1.1. Description

The WST-8 Cell Proliferation Kit is a colorimetric assay for the determination of viable cell number and for studying induction or inhibition of cell proliferation *in vitro*. This assay kit is based on the cellular reduction of the tetrazolium salt WST-8 [2-(2-methoxy-4-nitrophenyl)-3-(4-nitrophenyl)-5-(2,4-disulfophenyl)-2H-tetrazolium, monosodium salt] into a highly water-soluble, orange-colored formazan dye upon reduction in the presence of an electron carrier. As opposed to MTT assay, no solubilisation process is required since this formazan does not require solvation: the WST-8 is soluble in the tissue culture medium.

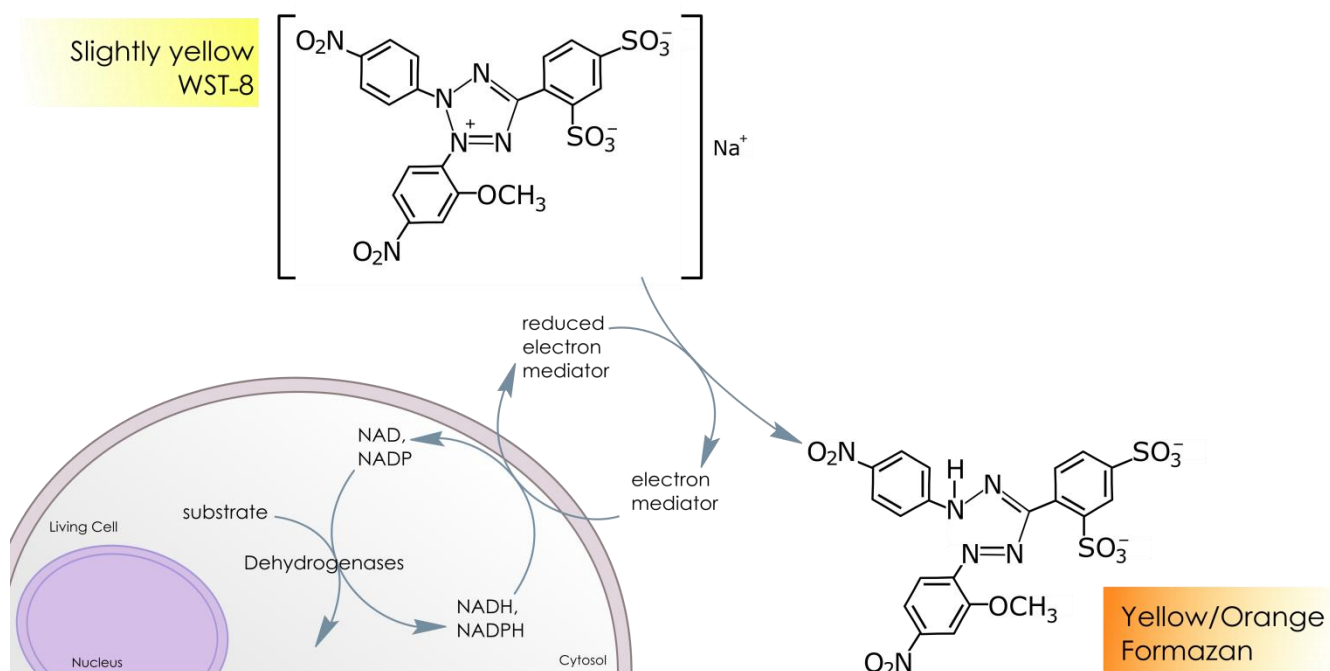

**Figure 1:** WST- 8 Cell Proliferation Kit mechanism: slightly yellow WST-8 is transformed into Yellow/Orange Formazan.

The WST-8 Cell Proliferation Kit is more sensitive at neutral pH than other tetrazolium salts such as MTT, XTT, MTS or WST-1 and the amount of formazan produced is directly proportional to the number of living cells.

Colorimetric measurement at 450 nm allows quantification of viable cells. The WST-8 Cell Proliferation Kit does not need cell fixation, cell lysis or washing steps rendering this ready-to-use kit fast, accurate, sensitive and adapted to high throughput screening.

## 1.2. Storage and shipping condition

**Storage:** Upon reception, store the WST-8 Cell Proliferation Kit at 4°C. Protect from light.

**Stability:** 1 year at 4°C

**Shipping condition:** The kit is shipped at RT.

## 2. Applications and Protocols

### 2.1. General Considerations

- Avoid repeated thawing and freezing cycles that could increase the background signal.
- This kit is compatible with phenol red containing medium.
- Incubation time varies with the type and number of cells; long incubation time (up to 4 hours) can be used for an optimal measurement.
- Allow reagent to reach room temperature before starting. Avoid direct exposure to- and protect from light.

### 2.2. General protocol for 96-well plate

#### *Cell preparation:*

1. Seed cells in a 96-well plate under standard culture conditions.
2. Carry out experiment by adding chemical compounds or biological agents to cells

#### *Measuring cell proliferation or viability:*

3. Add 10  $\mu$ L of WST-8 solution.
4. Incubate cells 30 min to 4h at 37°C in the dark (incubation time varies depending on the metabolic activity of the cells).
5. Optionally: centrifuge cells and apply protocol on the supernatant (see below)
6. Measure absorbance at 450 nm.
7. Subtract background absorbance of the non-treated cells from all other values.

Optionally for suspension cells, a centrifugation procedure to pellet cells (5 min x 1200 rpm) would allow working on supernatant to lower background signal.

### 2.3. General protocol for other well format


This kit is compatible with any cell culture format, from 384-well to 6-well plates. Follow the same protocol than for 96-well plates; changes only apply on the volumes used:

- Step 3: adjust the volume of the WST-8 solution in a well to 10% of the total volume.


## 2.4. Performance characteristics

### Cell titration and viability measurement after transfection.

This kit is compatible with both cell suspension and adherent cells. Cell titration was performed using JurKat T cell line cultivated in suspension and viability was measured using HEK-293 adherent cell line.

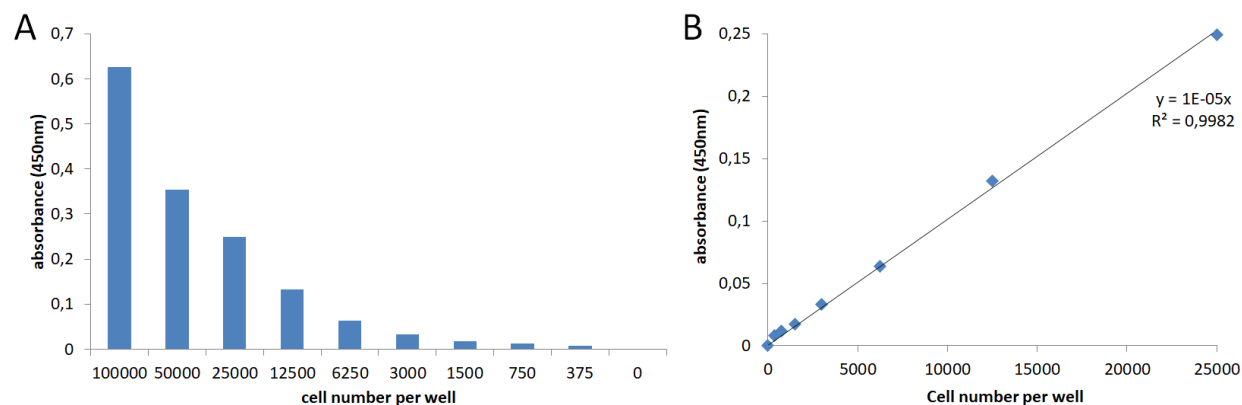

**Figure 2:** Cell number titration. (A) Jurkat T cells were seeded at different concentrations in a 96-well plate and incubated 2h with WST-8 Cell Proliferation Kit. Linearity between absorbance and cell number varies according to the cell type and metabolism (B).

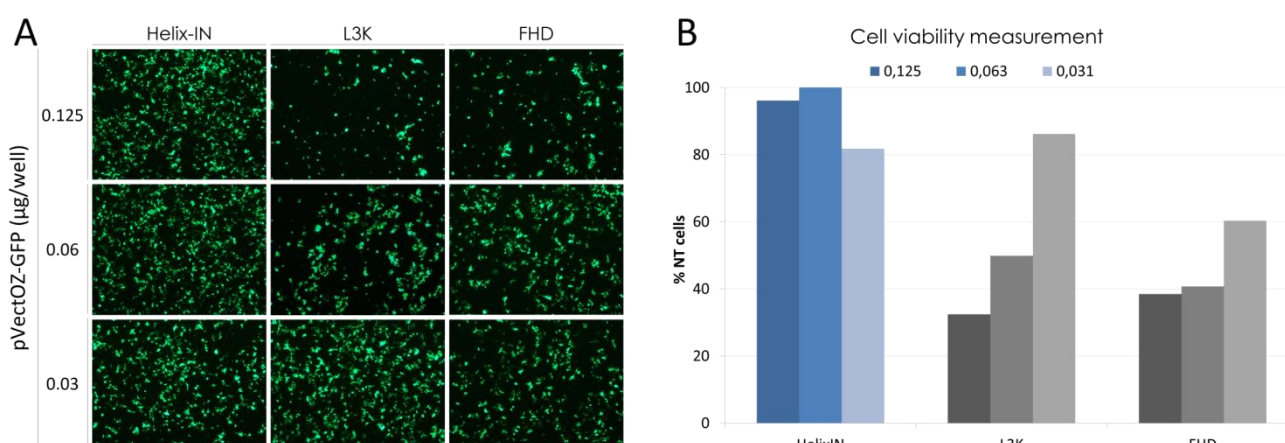

**Figure 3:** Cell viability measurement after transfection. (A) HEK-293 cells were cultivated in 96-well plates and transfected with 0.03, 0.06 and 0.125 µg DNA per well using Helix-IN DNA Transfection Reagent (OZ Biosciences) at a 2:1 ratio and two other transfection reagents according to their respective manufacturer's protocols. After 48h, transfection efficiency was monitored by fluorescence microscopy (B) and viability was determined as a % of Non-transfected (NT) cells using WST-8 Cell Proliferation Kit. 10µL of WST-8 Reagent were added to the cells and Absorbance at 450 nm was measured after 2h incubation at 37°C.

## **Purchaser Notification**

### **Limited License**

The purchase price paid for WST-8 Cell Proliferation Kit by end users grants them a non-transferable, non-exclusive license to use the kit and/or its separate and included components (as listed in the section 1, Kit Contents). These reagents are intended **for internal research only** by the buyer. Such use is limited to the use in the product manual. Furthermore, research only use means that this kit and all of its contents are excluded, without limitation, from resale, repackaging, or use for the making or selling of any commercial product or service without the written approval of OZ Biosciences.

Purchasers may terminate this License at any time by returning all WST-8 Cell Proliferation Kit material and documentation to OZ Biosciences, or by destroying all WST-8 Cell Proliferation Kit components. Purchasers are advised to contact OZ Biosciences with the notification that a WST-8 Cell Proliferation Kit is being returned in order to obtain a refund and/or to expressly terminate a research only license granted through the purchase of the kit(s).

This document covers in full the terms of the WST-8 Cell Proliferation Kit research only license, and does not grant any other express or implied license. The laws of the French Government shall govern the interpretation and enforcement of the terms of this License.

### **Product Use Limitations**

The WST-8 Cell Proliferation Kit and all its components are developed, designed, intended, and sold for research use only. They are not to be used for human diagnostic or included/used in any drug intended for human use. All care and attention should be exercised in the handling of the kit components by following appropriate research lab practices.

For more information, or for any comments on the terms and conditions of this License, please contact:

#### **Director of Business Development**

##### **OZ Biosciences SAS**

Parc Scientifique et Technologique de Luminy

Bâtiment Grand Luminy technopole

Case 922 zone entreprises

13288 Marseille Cedex 9, France

Ph: +33 (0)4.86.94.85.16

Fax: +33 (0)4.86.94.85.15

E-mail: [contact@ozbiosciences.com](mailto:contact@ozbiosciences.com)

## Related Products

PLASMIDS – REPORTER GENES
pVectOZ-LacZ pVectOZ-SEAP pVectOZ-GFP pVectOZ-Luciferase
ASSAY KITS
<b>Senescence / Viability / Stress</b> ROS Detection Assay Kit OZBlue Cell Viability Kit MTT Cell Proliferation Kit Senescence Kit for Stem Cells <b>Protein Dosage</b> FluoProdige Protein Quantification Assay Kit Bradford PaK BCA PaK <b>Reporter Gene</b> SEAP Assay Kit Luciferase Assay Kit $\beta$ -Galactosidase assay kits (CPRG/ONPG/X-Gal Staining Kit)

## CONTACTS

**OZ Biosciences SAS**  
163 avenue de Luminy  
Case 922, zone entreprise  
13288 Marseille cedex 09  
**FRANCE**

Ph: +33 (0) 486 948 516  
Fax: +33 (0) 486 948 515

[contact@ozbiosciences.com](mailto:contact@ozbiosciences.com)  
[order@ozbiosciences.com](mailto:order@ozbiosciences.com)  
[tech@ozbiosciences.com](mailto:tech@ozbiosciences.com)

**OZ Biosciences INC**  
4901 Morena Blvd,  
Suite 501  
San Diego CA 92117  
**USA**

Ph : + 1-858-246-7840  
Fax : + 1-855-631-0626

[contactUSA@ozbiosciences.com](mailto:contactUSA@ozbiosciences.com)  
[orderUSA@ozbiosciences.com](mailto:orderUSA@ozbiosciences.com)  
[techUSA@ozbiosciences.com](mailto:techUSA@ozbiosciences.com)

<http://www.ozbiosciences.com>

Follow us!

