



< Kaneka Easy DNA Extraction Kit version 2* >

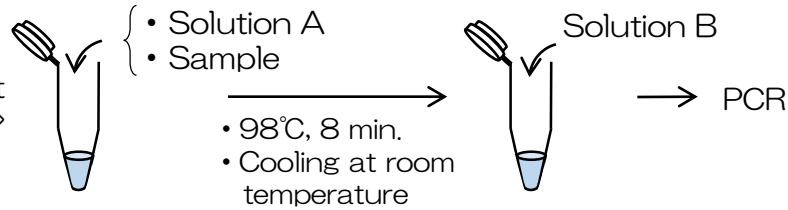

Simple protocol without centrifugation

< A conventional method >

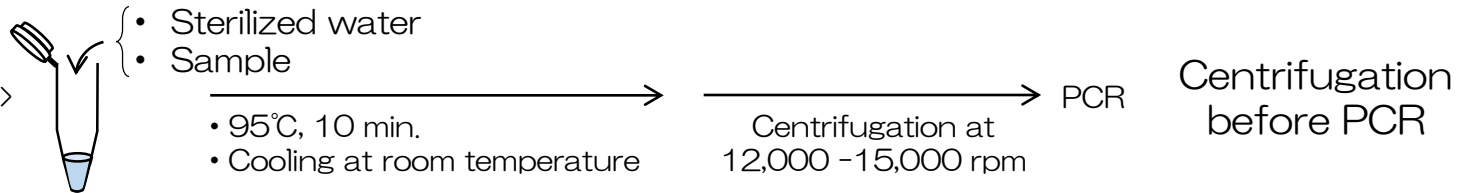

< DNA extraction from candida colony and following PCR >


1. Pick up candida colonies in solid media with inoculation loops.
2. Extract DNA from candida colonies by using Kaneka Easy DNA Extraction Kit version 2.



PCR and electrophoresis for confirming DNA amplification

(Data obtained by our company)

M: 50 bp DNA ladder marker
 N: negative control (water)

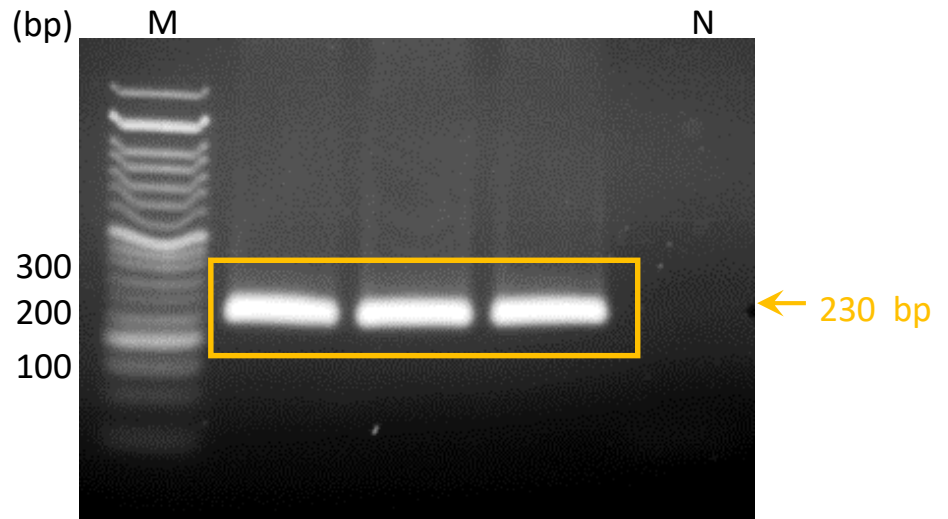

Kaneka DNA Easy Extraction Kit version 2 successfully extracted DNA from candida colonies.

* "Kaneka Easy DNA Extraction Kit version" 2 is for research use only.