

Kaneka Easy DNA Extraction Kit Version 2 [Research Use Only]

For more information : http://www.funakoshi.co.jp/exports_contents/80099

Easy DNA extraction kit for routine template DNA

2 reagents and rapid around 10 minutes protocol

Extraction kit for cell lysates, animal and plant tissues, seed, bacteria, blood and feces



Features

- Simple protocol
- Time saving - Around 10 minutes per sample for extraction
- Room Temperature Storage
- Broad range of sample types are compatible with this kit
 - Cell lysates, Animal tissues, Plant tissues, Seed, Bacteria, Blood and Feces
- 1 kit includes 250 preps (for Blood sample) or 50 preps (for other samples) of reagents


Example Protocol

(For Plant Seeds Sample)

1. Put a 5 ~ 8 mm square piece of a seed (cut by scissors etc.) into a PCR tube. Add 100 μ L of Solution A to the tube and mix well by pipetting.
2. Incubate the PCR tube at 98°C for 8 minutes in a heat block or other heating equipment.
3. After the PCR tube is cooled, add 14 μ L of Solution B and mix well.
4. Stir the extract from Protocol 3 above well before use. Use 1 ~ 5 μ L of it as a template DNA for PCR (for a 50 μ L PCR reaction).

Protocol for other samples are also available !


Results

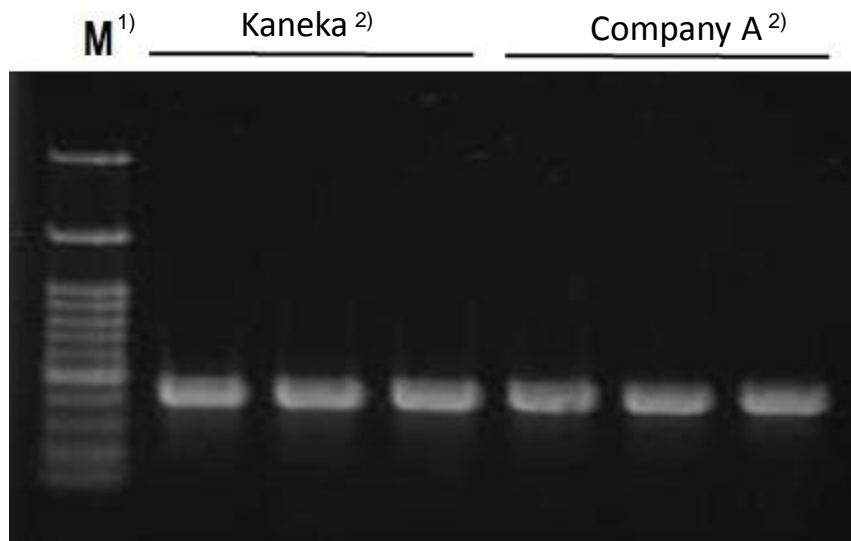

1) DNA ladder marker
2) Number of sample: n=3

Extraction from mouse tail sample

DNA were extracted from mouse tail sample with Kaneka kit or Company A kit and extracted DNA samples were analyzed by end point PCR.
PCR data showed the same efficiency in extraction as Company A kit (around 13 minutes protocol).

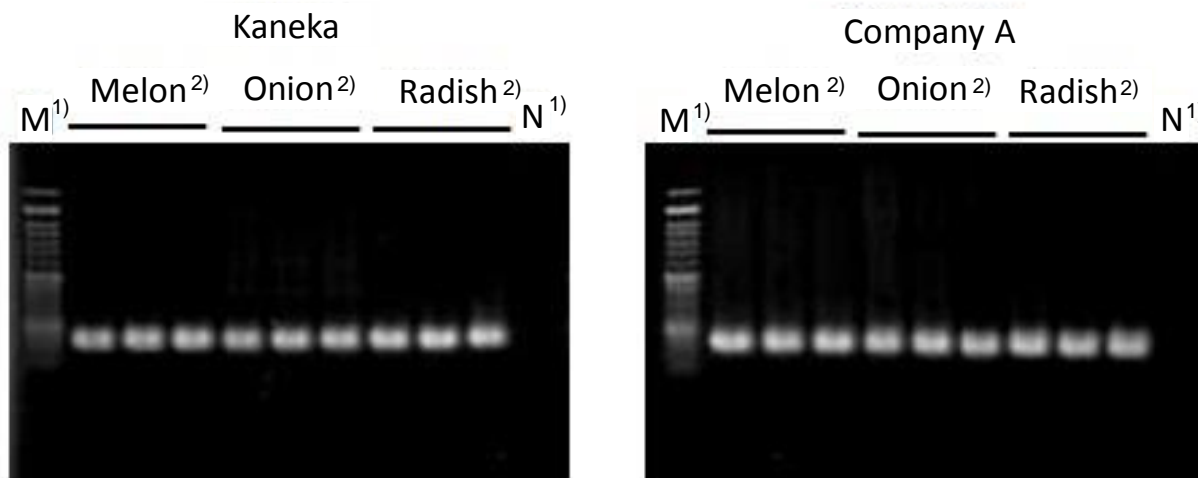

1) DNA ladder marker 2) Number of sample: n=3

Extraction from plant seed samples

DNA were extracted from plant seed samples (Melon, Onion, or Radish) with Kaneka kit or Company A kit and extracted DNA samples were analyzed by end point PCR.
PCR data showed the same efficiency in extraction Company A kit (around 80 minutes protocol).

Product Information

[Manufacturer : KNK]

Product Name	Size	Catalog #	Storage
Kaneka Easy DNA Extraction Kit Version2	1 kit	KN-T110005	RT

NOTE

※ All products here are research use only, not for diagnostic use.
※ Specs might be changed for improvement without notice.

※ Company name and product name are trademark or registered mark.
※ Please contact your local distributors for orders, quote request and inquiry.

Your Local Distributor

funakoshi Co., Ltd.

Address: 9-7 Hongo 2-Chome, Bunkyo-ku,
Tokyo 113-0033 JAPAN
Phone : +81-3-5684-6296
Fax : +81-3-5684-6297
Email : export@funakoshi.co.jp

ZF-Y07T-08 (2015.08)