



Premier BOX-V19 Example of use

《Set material》

- iP-TEC Temperature Stabilizer-36 6pcs
 - Secondary Container of Live-Transport 2set
 - iP-TEC Flask-25 12pcs
- Application using the above

- iP-TEC Temperature Stabilizer-36 8pcs
 - Secondary Container of Live-Transport 4set
 - iP-TEC Flask-25 24pcs
- Application using the above



Spacer
2pcs



Band Cushion
4pcs



Mesh Cushion
4pcs



Set two spacers and four Band Cushions as shown in the photograph.



Place two Temperature Stabilizer on the bottom and set.




Put the first set of secondary containers. Also set Temperature Stabilizer on the side (put on the band cushion). Fix the secondary container with 2 mesh cushions.




Then put one Temperature Stabilizer on top of it.




Place the second secondary container set on top of it. Then fix the secondary container with two mesh cushions.

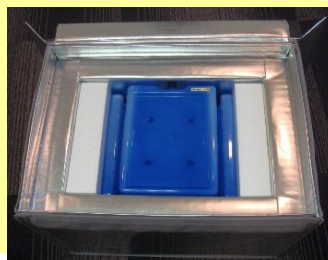

Place a piece of Temperature Stabilizer and cover it.

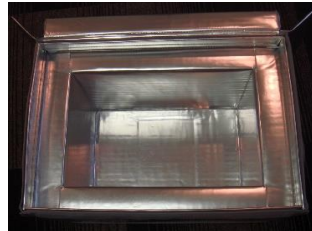


With two Temperature Stabilizer stacked, set two columns.

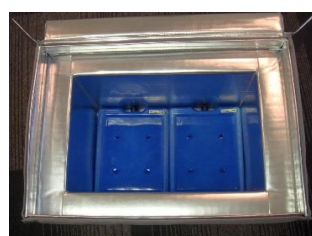

Place the secondary container set on top of each. Fix the secondary container with 4 band cushions.



One Temperature Stabilizer is placed on each

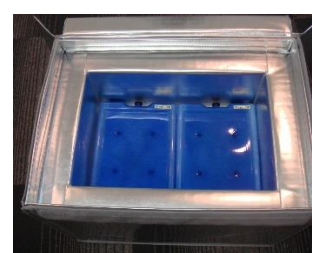

Place the third and fourth secondary container set on top of it. Fix the secondary container with 4 mesh cushions.



Place Temperature Stabilizer on each and cover it.

* Storage of premium BOX when unused is best under low temperature / low humidity environment.



This product is for research use, it is not a specification assuming medical use.



SANPLATEC CORP.